


# TEMVu™

Electron Microscopy Meets Light Microscopy

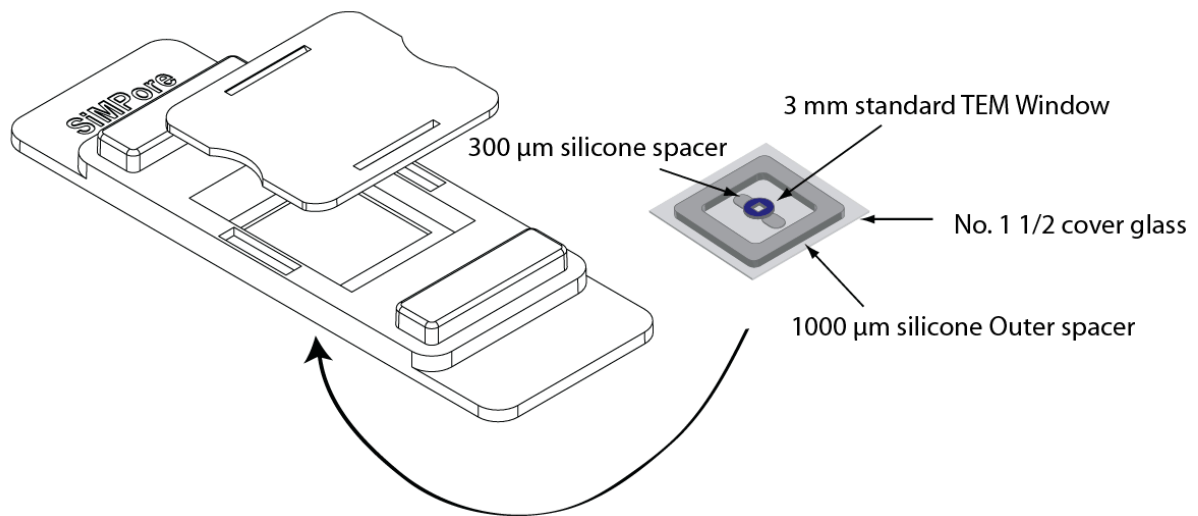

Ultrathin Membranes for Light, Fluorescent and  
Electron Microscopy

TEMVu™ is for research use only  
and not for use in diagnostic or clinical applications.

**Customer Service**

www.TEMwindows.com  
service@temwindows.com  
phone: 585-214-0585

**Order Information**

www.TEMwindows.com  
sales@temwindows.com  
phone: 888-323-NANO  
fax: 888-249-2935



## OPERATIONS

### Required Equipment

- ◆ Pipette (2-200  $\mu$ L tip)
- ◆ Autoclave for sterilization (for sterility)
- ◆ Standard cell culture media and incubation equipment (for cell culture on TEMVu™)
- ◆ Standard fixatives/preferred stains.

### Instructions

Please read all instructions and notes before proceeding

#### Cell Culture Instructions

1. Remove slide carrier from plastic bags and remove inspection label. Do NOT press glass window. The glass is thin for imaging purposes and therefore delicate.
2. Sterilize TEMVu slide and cap by autoclaving for cell culture applications
3. To improve cell adhesion, incubate TEMVu grid with 10  $\mu$ L cell media for at least two hours prior to introducing cells. Gently aspirate off this incubation media just prior to seeding cells. Membranes can also be exposed to ECM proteins in solution prior to cell seeding.
4. To wet the bottom of the TEMwindow grid, pipette  $\sim$ 10  $\mu$ L of cell culture fluid using a fine pipette tip, under the TEMwindow grid. This will enable both sides of the membrane to stay wet and will help cell growth and imaging quality.
5. To seed cells on the TEMwindow grids, pipette 10  $\mu$ L of suspended cells ( $4\text{-}8 \times 10^5$  cells/mL) onto the top of TEMwindow grid to allow cells to settle on to the membrane surface.
6. After two (or more) hours of incubation, the cells should have attached to the membrane surface. Gently aspirate off media to remove any unattached cells. Load fresh culture media ( $\sim$ 150 $\mu$ L) into the chamber.
7. Replace the TEMVu cap and place directly in the incubator or within a petri dish. TEMVu is designed to be stackable to preserve space.

#### Staining/Imaging Instructions

1. Cells/Materials deposited on the grids can be visualized at any point using upright or inverted microscopes. The grids do not show autofluorescence and will not interfere with imaging.
2. Staining and fixing of cells can be performed directly on TEMVu slide if needed. The silicone gasket has limited compatibility with organic solvents.
3. Once staining/fixing is complete, the TEMwindow grids can be removed using the included K6 plastic tweezers (cat no. K6TWZR).

#### NOTE:

- Due to the extreme thinness of TEMVu™ membranes it is important to aspirate as gently as possible. Do not directly touch the membrane with a pipette tip. Vacuum aspiration may be too harsh in some laboratory environments. In these cases, gently use a hand pipette.

- If you have trouble wetting the underside of the TEMwindow grid, tilt the slide away from the pipette tip before introducing the fluid. A fine tip pipette such as gel loading tip can also be used. If this approach is not working, very gently lift the grid with the plastic tweezers to help wet the underside.
- It is important to keep the both sides of membranes wet for cell culture, as cells tend to grow differently if one side of the membrane is dry.
- To achieve optimum image quality, make sure underside of TEMwindow grid is wet and free from air bubbles.
- If silicone gasket/plastic incompatible solvents have to be used, the TEMwindow grids can be gently removed from the TEMVu slide using the included K6 plastic tweezers and processed further.

If you have any additional question please do not hesitate to contact us.

## SPECIFICATIONS

### Sterilization

TEMVu™ slides can be sterilized by autoclaving after removing from plastic bag and discarding the inspection label.

### Chemical Stability

Devices are incompatible with strong bases. Silicon based nanoporous membranes degrade over a period of 3 days in physiological buffers (pH 7 or greater). Non-degradable nanoporous membranes (available upon request) are stable for at least 2 weeks, while nitride based membranes are stable for at least 4 weeks in physiological buffers. The dissolution of the membrane is non-toxic and has not been shown to adversely affect cell viability. If using solvents, please verify chemical compatibility against silicone gasket and polycarbonate slide.

### Device Storage

Store in a clean and dry environment. Exposure to UV can damage polycarbonate slide carrier.


### TEM Grid Specifications

TEMVu™ Slides are built to order. All standard TEMwindow grids available at [www.TEMwindows.com](http://www.TEMwindows.com) can be assembled into your slides. Please visit the website for a complete listing of grids available. Slides will ship within 5 days of your order.


### Physical Properties

<b>Product Number</b>	TEMVU-(Individual Grid Product Number)
<b>Membrane Compositions Available:</b>	Porous Nanocrystalline Silicon (pnc-Si), Non Porous Silicon, Silicon Dioxide or Silicon Nitride
<b>Stability in Culture:</b>	3 days for Porous Nanocrystalline Silicon (pnc-Si), Non Porous Silicon, Silicon Dioxide ~ 2 weeks for Non-degradable Porous Silicon ~4 weeks for Silicon Nitride
<b>Dimension of TEM Grid:</b>	3 mm Standard Window
<b>Chamber Volume:</b>	150 µL
<b>Glass Support:</b>	No. 1 ½
<b>Slide Composition:</b>	Polycarbonate
<b>Gasket Composition:</b>	Medical Grade Silicone
<b>Gasket Thickness:</b>	100/300 µm

## SCHEMATICS

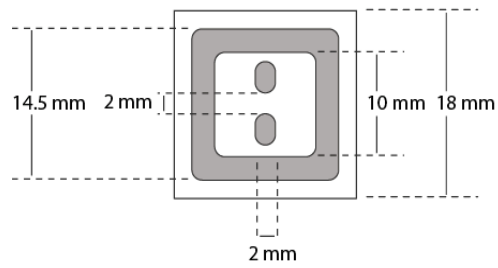

## Cross Sectional View

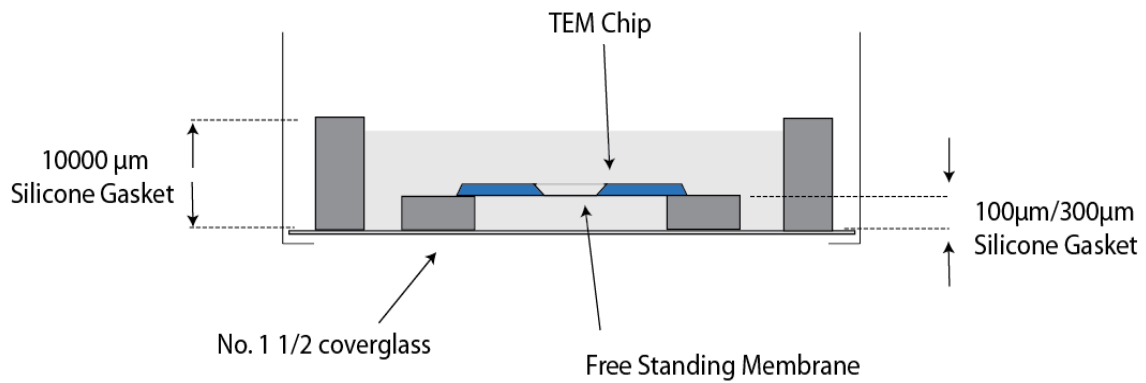

## STANDARD WARRANTY

**SiMPore Inc.** ("SiMPore") warrants its products will meet their applicable published specifications when used in accordance with their applicable instructions for a period of one year from shipment of the products. **SIMPORE MAKES NO OTHER WARRANTY, EXPRESSED OR IMPLIED. THERE IS NO WARRANTY OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE.** The warranty provided herein and the data, specifications, and descriptions of SiMPore products appearing in SiMPore's published catalogues and product literature may not be altered except by express written agreement signed by an officer of SiMPore. Representations, oral or written, which are inconsistent with this warranty or such publications are not authorized and if given, should not be relied upon.

In the event of a breach of the foregoing warranty, SiMPore's sole obligation shall be to repair or replace, at its option, the applicable product or part thereof, provided the customer notifies SiMPore promptly of any such breach. If after exercising reasonable efforts, SiMPore is unable to repair or replace the product or part, then SiMPore shall refund to the customer all monies paid for such applicable product or part. **SIMPORE SHALL NOT BE LIABLE FOR CONSEQUENTIAL, INCIDENTAL, SPECIAL, OR ANY OTHER INDIRECT DAMAGES RESULTING FROM ECONOMIC LOSS OR PROPERTY DAMAGE SUSTAINED BY ANY CUSTOMER FROM THE USE OF ITS PRODUCTS.**

Version 1106